

Compact Photodiode Power Head with Silicon Detector



S121C

Description

The S121C power head is designed for general purpose optical power measurements. The head is optimized for small thickness to fit in tight spaces. The high sensitive photodiode with large active area in combination with a reflective, diffused ND filter enables power measurements up to 500 mW in free-space and fiber-based applications. A removeable annular VIS/IR viewing target allows conveniently centering the measured beam to the active area of the photo-diode. The target absorbs light from 400 to 640nm and 800 to 1700nm.

The S121C housing includes a threaded input in axis with the light input aperture that is compatible with any number of Thorlabs 1" threaded accessories. This allows convenient mounting of external optics, fiber adapters, light filters, and apertures. A #8-32 and M4 threaded mounting hole is provided to accommodate posts and post holders.

The S121C is compatible with all new Thorlabs display units. A non-volatile memory in the sensor connector contains sensor information data and the NIST and PTB traceable calibration data.

Available Accessories

| | |
|----------|---------------------|
| S120-FC | FC fiber adapter |
| S120-SMA | SMA fiber adapter |
| S120-SC | SC fiber adapter |
| S120-LC | LC fiber adapter |
| S120-ST | ST fiber adapter |
| SM1CP1 | Dust cap (included) |

The S121C is also compatible to the Thorlabs imperial and metric post and post-holder series and Thorlabs SM1 mechanics.

Cleaning and Maintenance

There are no serviceable parts in the S121C head. The housing may be cleaned by wiping with a soft damp cloth. When cleaning the aperture filter, treat it as any other fine optic. Gently blow off any debris using compressed air and wipe gently with an optic tissue wetted with propanol. If you suspect a problem with your S121C please call Thorlabs and an engineer will be happy to assist you.

As long as the sensor has not been exposed to excessive optical power (please pay attention to the maximum ratings in the technical specifications), the calibration should be very stable over long periods of time (well over a year). To keep the accuracy and performance of the S121C, Thorlabs recommends a yearly recalibration, starting one year after purchase.

Specifications

| | |
|---------------------------------------|---|
| Detector Type | Silicon Photodiode |
| Wavelength Range | 400 - 1100 nm |
| Optical Power Working Range | 500 nW - 500 mW |
| Max Average Power Density | 20 W/cm ² |
| Max Pulse Energy | 20 μJ |
| Linearity | ± 0.5% |
| Resolution ¹⁾ | 10nW |
| Measurement Uncertainty ²⁾ | ±3% 440 - 980 nm ±5% 400 - 439 nm, ±7% 981 - 1100 nm |
| Typical Application | Low Power Lasers |
| Laser Types | Diode, Diode Arrays, He-Ne, Dye, Ion Lasers (Ar+, Kr+) |
| Coating /Diffuser ³⁾ | Reflective ND (OD2) |
| Cooling | Convection |
| Head Temperature Measurement | NTC Thermistor 4.7kΩ |
| Console Compatibility | PM100D, PM100A, PM100USB, PM200, PM320E |
| Response Time | < 1 μs |
| Sensor Dimensions | Ø30.5 mm x 12.7 mm |
| Active Detector Area | 9.7 mm x 9.7 mm |
| Input Aperture | Ø9.5 mm |
| Cable Length | 1.5 m |
| Connector | Sub-D 9p male |
| Weight | 0.07 kg |
| Post ³⁾ | #8-32 & M4 thread |
| Aperture Thread | SM1, outer thread |
| Fiber Adapters (optional) | FC, SC, LC, SMA, ST |


¹⁾ Measured with PM100D console in bandwidth low setting.

²⁾ Beam diameter > 1mm


³⁾ This specification is valid for S121C devices from serial number 1203xxx. For older versions, please contact technical support.

Please note that the S121C power meter head is not compatible with the older Thorlabs power meter consoles (PM100, PM30, PM300, PM300E, S100).

Typical Response Graph

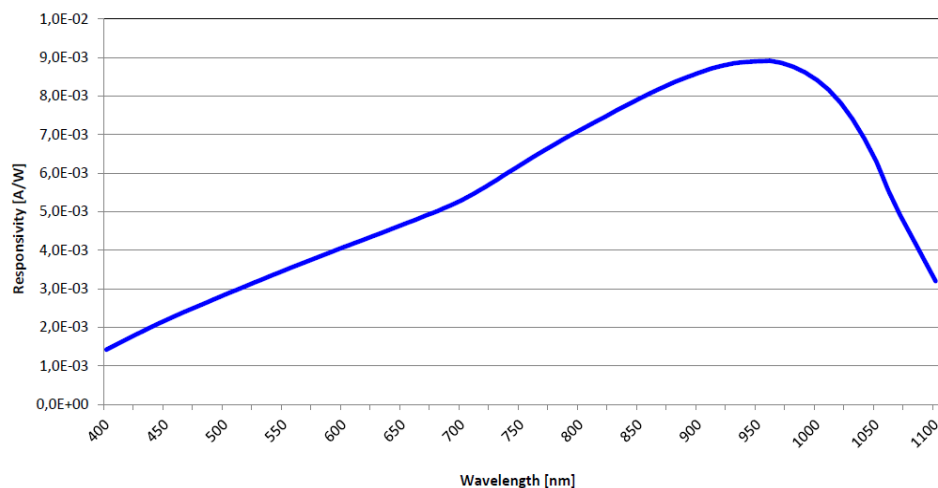

Drawings

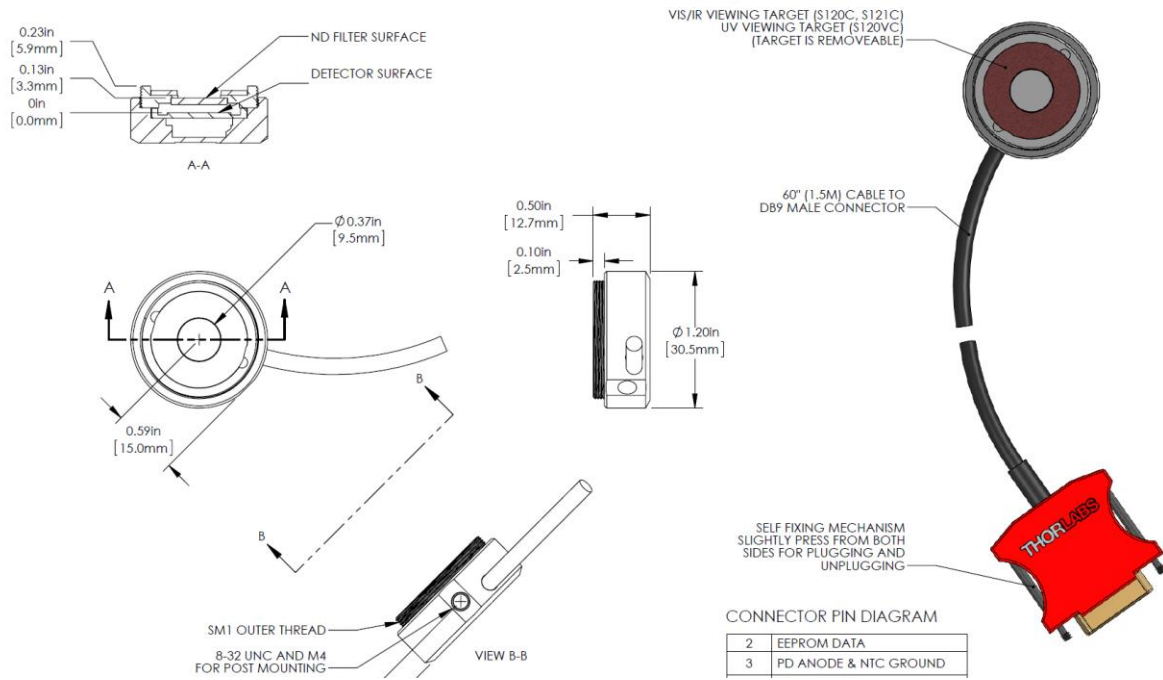

Precautions and Warranty Information

These products are ESD (electro static discharge) sensitive and as a result are not covered under warranty. In order to ensure the proper functioning of a photodiode care must be given to maintain the highest standards of compliance to the maximum electrical specifications when handling such devices. The photodiodes are particularly sensitive to any value that exceeds the absolute maximum ratings of the product. Any applied voltage in excess of the maximum specification will cause damage and possible complete failure to the product. The user must use handling procedures that prevent any electro static discharges or other voltage surges when handling or using these devices.

Thorlabs, Inc. Life Support and Military Use Application Policy is stated below:

THORLABS' PRODUCTS ARE NOT AUTHORIZED FOR USE AS CRITICAL COMPONENTS IN LIFE SUPPORT DEVICES OR SYSTEMS OR IN ANY MILITARY APPLICATION WITHOUT THE EXPRESS WRITTEN APPROVAL OF THE PRESIDENT OF THORLABS, INC. As used herein:

- 1. Life support devices or systems are devices or systems which, (a) are intended for surgical implant into the body, or (b) support or sustain life, and whose failure to perform, when properly used in accordance with instructions for use provided in the labeling, can be reasonably expected to result in a significant injury to the user.*
- 2. A critical component is any component in a life support device or system whose failure to perform can be reasonably expected to cause the failure of the life support device or system or to affect its safety or effectiveness.*
- 3. The Thorlabs products described in this document are not intended nor warranted for usage in Military Applications.*

